# Fagron Sterile (FAST)

**The Aseptic Solution**

# Introduction

Ger van Jeveren  
Founder and CEO

UK

## The FAST Aseptic Solution

Fagron provides over 200,000 customers worldwide with high-quality concepts and innovations, making it possible to provide tailor-made pharmaceutical care to patients.

Our ability to listen to our customers and work closely with prescribers, pharmacists and academics is the foundation of our successful strategy of optimising and innovating pharmaceutical compounding.

With a continuous and fast growing pipeline of innovations, 150 pharmacists dedicated to compounding and a range of professional training programmes, Fagron widens the therapeutic scope of the Prescriber and Compounding Pharmacist to offer bespoke pharmaceutical solutions.

Now... Fagron has worked closely with specialists and suppliers to develop and introduce a new concept that incorporates an extensive range of equipment and materials for the Aseptic Compounder.

### We call this our FAST concept.

The FAST concept uniquely combines leading-edge equipment with minimum validation requirements and innovative accessories to help reduce time-consuming repetitive manual operations and the opportunity for human error.

Developed to save time and laboratory space our easy to operate, minimum footprint single and multiple channel closed system pumps accurately deliver mixed infusion solutions on time... every time.

Combined with our revolutionary Fagron Aseptic Pack (FAP) infusion bags and an unrivalled portfolio of easily available essential accessories and consumables, the FAST concept is without competition.

### Fast, Accurate, Sterile and on Time.

Making life easier and more efficient for the Compounder and ultimately much better and safer for the Patient.

The comprehensive range of products in our FAST service can be found in the Fagron FAST catalogue.

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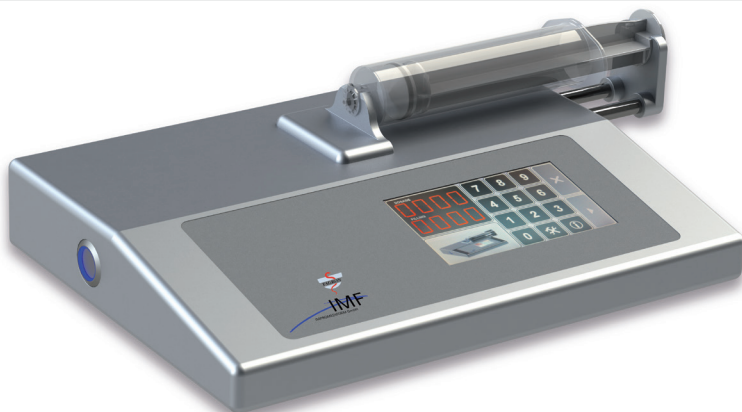
UK

## Features/Benefits:

- Low installation and validation costs
- Lightweight, easy to move
- Low noise emission
- Simple to operate and easy to clean
- High speed filling - up to 220ml max./per minute
- Selectable programmes - e.g. for sterile filtration
- 12 month warranty (extendable), with option for service contract
- Single or multiple syringe filling possible
- For mixing, transferring and filling liquids quickly and accurately.

For example:

- Total parenteral nutrition (TPN)
- Filling devices (including vials, syringes and elastomeric pumps)
- Chemotherapy, antibiotics
- Patient-specific infusions
- Filling syringes
- Paediatric doses
- Sterile filtration

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Single channel volumetric pump for mixed infusion solutions in a closed system delivery.

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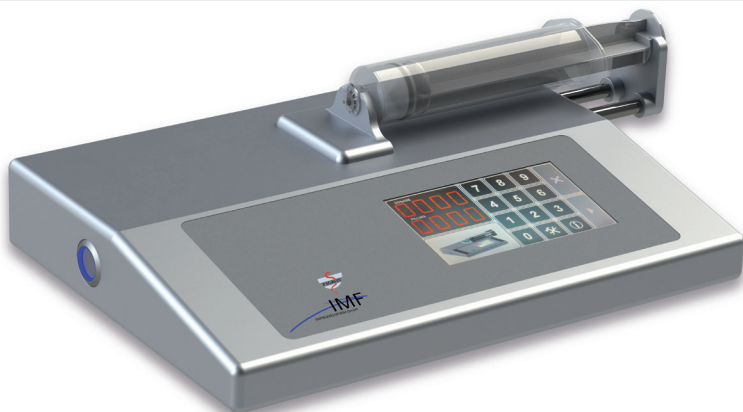
UK

## Technical Data:

Syringe volume:	50ml
Max pumping rate:	220ml/min
Input accuracy:	1ml steps; optionally 0.1ml
Volume entry range:	5 - 9999ml
Accuracy:	metering volumes greater than 10ml: less than 1%; metering volumes 5 to 10ml: less than 3%
Power supply:	100-240V, 50/60Hz
Power:	Max. 25VA
Size (WxDxH):	280 x 205 x 50mm
Space required (WxDxH): (including filled syringe)	400 x 205 x 60mm
Weight:	3.5kg
Noise emission:	< 60dB(A)

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## MF4014

UK

### Paediatric Metering System (4 Supply Lines)

4 ways-in system for the 1-Channel Pump

For precise millilitre admixturing of TPN, infusion therapy and liquids in a closed system

- valve operated
- easy handling
- 4 lines
- can be increased to 7 lines
- bacteria-dense aeration
- integrated 15µm particle filters
- drip chamber as a solution reservoir
- 50ml dosage syringe
- different coloured large tubing clamps
- rotatable luerlock thread nut

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UK

## Features/Benefits:

- Low installation and validation costs
  - Lightweight, easy to move
  - Low noise emission
  - Simple to operate and easy to clean (gassable)
  - High speed filling - up to 220ml max./per minute per channel
  - Selectable programmes - e.g. for sterile filtration
  - Simple to operate with integrated remote tablet control screen
  - Barcode reader reduces risk
  - Optional link to formulation software
  - 12 month warranty
  - Single or multiple syringe filling possible
  - For mixing, transferring and filling liquids quickly and accurately
- For example:
- Total parenteral nutrition (TPN)
  - Patient-specific infusions
  - Paediatric doses
  - Chemotherapy, antibiotics
  - Filling devices (including vials, syringes and elastomeric pumps)
  - Filling syringes
  - Sterile filtration

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Six channel volumetric pump for multi-source fluid mixing in a closed system delivery.

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UK

## Technical Data:

<b>Syringe volume:</b>	50ml
<b>Max pumping rate:</b>	220ml/min per channel
<b>Input accuracy:</b>	1ml steps; optionally 0.1ml
<b>Volume entry range:</b>	5 - 9999ml
<b>Accuracy:</b>	metering volumes greater than 10ml: less than 1%; metering volumes 5 to 10ml: less than 3%
<b>Power supply:</b>	100-240V, 50/60Hz
<b>Power:</b>	Max. 150VA
<b>Size (WxDxH):</b>	520 x 240x 400mm
<b>Space required (WxDxH):</b>	275 x 220 x 45mm
<b>(including filled syringe without tablet interface)</b>	
<b>Weight:</b>	13kg/2.6kg
<b>Noise emission:</b>	< 65dB(A)

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## UK EVA Infusion Bags with Luerlock

Lf St Nd

Sterile EVA bag with luerlock connector. Non Latex additive port. Add port, filling port and giving port.

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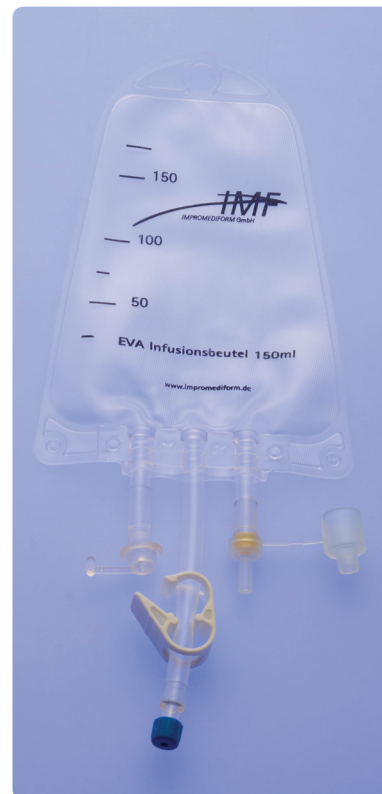
## UK EVA Infusion Bags with Luerlock - Light Protected

Lf St Nd

Sterile EVA bag with luerlock connector. Non Latex additive port. Add port, filling port and giving port. Light protected.

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## UK EVA Infusion Bags - Big Bore

Lf St Nd

Sterile EVA bag with large-bore threaded connector. Non Latex additive port. Add port, filling port and giving port.

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## UK EVA Infusion Bags - Light Protected Big Bore

Lf St Nd

Sterile EVA bag with large-bore threaded connector. Non Latex additive port. Add port, filling port and giving port. Light protected.

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# EVA Total Parenteral Nutrition Bags

## UK **EVA Infusion bags with 3-way Transfer Sets**

Lf St Nd

Sterile EVA bag with large-bore threaded connector. Non Latex additive port. Add port, filling port and giving port. 3 Vented transfer spikes for gravity filling.

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## UK **EVA Dual-Chamber Bags for TPN**

Lf St Nd

Sterile EVA dual Chamber bag with luerlock connectors. Non Latex additive port. Add port, filling port and giving port.

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UK

### 3-Way Feeding Set

For transferring liquids from containers (e.g. glass bottles) by gravity into Infusion Bags MF1661 to MF1666

- 3 spikes
- large tubes
- high volume
- easy handling
- bacteria-dense aeration
- different coloured tubing clamps
- rotatable luerlock thread nut

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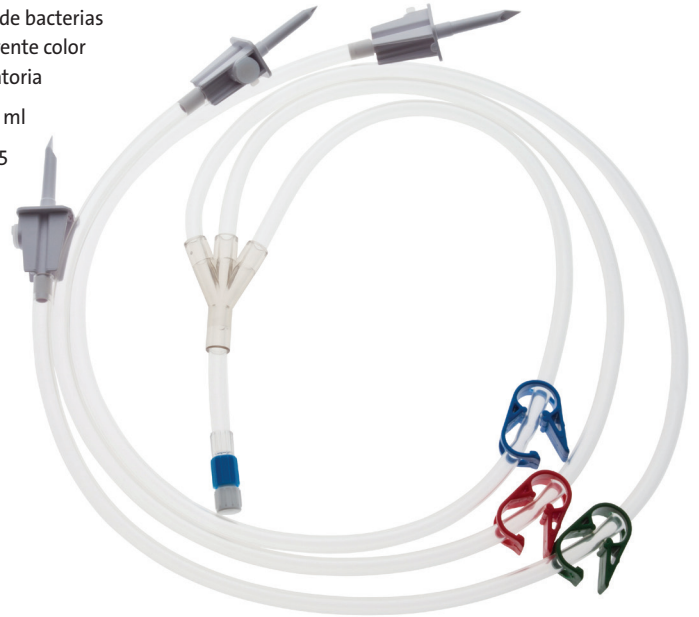
## MF8881

UK

### Infusion Line for Taxol® with 0.2µm Bacteria Filter

- PVC-free
- tube is made of MPE
- 0.2µm Supor infusion filter

The infusion line is pressure-proof and free of PVC. The recommended 0.2µm filter for administering Taxol consists of an air separator and is provided with a 0.2µm Supor membrane

The infusion tube is made of Polyethylene (Metalloccen Polyethylene: MPE) which does not keep back the Paclitaxel agent. The retaining value is practically negligible

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UK

### Mixing Adaptor Flat

For the mixing and transfer of substances from Vials into Viaflex® bags

- large cannula
- minimal risk of blockage
- fixed needle with tapered needle holder
- increased patient and user safety

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### Mini Spike

SmartSite® needle-free system

Vented aspiration spike for liquid extraction from storage bottles and vials of various sizes.

- anti-aerosol
- 0.2µm hydrophobic microporous GORE-TEX®
- single packed

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### Twin Spike Valve

SmartSite® needle-free system

Vented aspiration spike for liquid extraction from storage bottles and vials of various sizes

- anti-aerosol
- 0.2µm hydrophobic microporous GORE-TEX®
- single packed

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UK

### Vented Spike

For puncturing stoppers and membranes, withdrawing or filling glass bottles or other containers

- bacteria-dense aeration (can be closed for infusion bags)
- tubing clamp

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# Fagron Aseptic Pack

## UK Efficient, easy and safe...

Preparing aseptic medication is a complex process. During the preparation of parenteral medication critical steps need to be taken. The opportunities for preparation errors are high. High risks such as microbiological contamination and patient harm due to mistakes in preparation are examples of these possible risks.

### How safe and fast is your aseptic preparation?

### Fagron offers you the right solution! Meet FAP...

Fagron Aseptic Pack is a unique concept - an easy and safe way to compound a tailor-made aseptic preparation. The weight and quantity of both API and diluent are guaranteed and, when combined in the dual-compartment Fagron Aseptic Pack contribute to optimal medication safety. The aseptic preparation is ready for use in just 3 steps.

### Fagron Aseptic Pack is efficient, easy and safe.

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**1: Presser, 2: Mélanger, 3: Préparer ...**

## UK Step 1: Press

Place the FAP flat on a firm surface, press with both hands on the diluent compartment to break the seal. Alternatively, using one corner of the diluent compartment gently roll the bag applying pressure on the seal until the seal breaks.

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# Fagron Aseptic Pack

## UK Step 2: Mix...

Mix the contents by inverting the FAP several times and ensure the API has completely dissolved.

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## UK Step 3: Compound...

Use a MediMix pump and the Vented Spike (MF1512) to fill the administration device (for example a syringe), or if filling small quantities by hand, attach to the FAP to compound the finished product.

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UK

## FAP advantages

- Saves resources - time and money and manpower
- Easy to use
- Enhances medication safety
- Safe and hygienic to work with
- Reduces/removes Repetitive Strain Injuries (RSI) risk
- Frees up the cleanroom for other activities
- Easy to store, wipe, clean and transport
- Long shelf life
- Efficient, easy and safe!

## FAP facts

- Weighed quantities of API and diluent in separate compartments
- Universal connection possibilities, including automated conditioning equipment
- Double vial stoppers for maximum flexibility for the tailor-made preparation
- No spillage due to unique design
- Latex-free and PVC-free

## FAP Individual Tailor-made Care

To provide optimal pharmaceutical care, tailor-made medication is of great importance. Fagron supports customers with individual tailor-made care - with FAP.

FAP individual tailor-made care offers solutions for specific needs from the pharmaceutical market, and Fagron ensure that FAP innovations fit in perfectly with our customers' specific requirements.

**If you have a product suggestion for a tailor-made FAP to make life easier for you, contact us and we will be happy to discuss your requirements.**

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# Index

UK

Product Code	Description	Page	Product Code	Description	Page
MF1504	Double Luerlock (Female)	11	MF1866	EVA Infusion Bag 3000ml with 3-way Transfer Set	22
MF1512	Vented Spike	35	MF1873	3-Way Feeding Set	23
MF1516	Transfer Set	10	MF4010	MediMix mini Pump	5
MF1544	SmartSite®	31	MF4011	Transfer System 1-Way	7
MF1590	Mixing Adaptor Flat	32	MF4012	Dosage System with 2 Lines	8
MF1593	Mini Spike	33	MF4014	Paediatric Metering System (4 Supply Lines)	9
MF1594	Twin Spike Valve	34	MF4060S	MediMix plus Pump	15
MF1605	Connection Pipe 15cm	12	MF4061/4062	Dosage System with 6 Lines into 6 Syringes (Horizontal/Vertical)	18
MF1606	Connection Pipe 30cm	13	MF4064	Dosage System with 4 Lines into 4 Syringes	17
MF1629	Syringe Pump Line 250cm	14	MFL516	Transfer Set Light Protected	10
MF1660	EVA Infusion Bag 100ml with Luerlock	20	MFL660	EVA Infusion Bag 100ml with Luerlock Light Protected	20
MF1661	EVA Infusion Bag 150ml with Luerlock	20	MFL661	EVA Infusion Bag 150ml with Luerlock Light Protected	20
MF1661B	EVA Infusion Bag 150ml Bigbore	21	MFL661B	EVA Infusion Bag 150ml Light Protected Bigbore	21
MF1662	EVA Infusion Bag 250ml with Luerlock	20	MFL662	EVA Infusion Bag 250ml with Luerlock Light Protected	20
MF1662B	EVA Infusion Bag 250ml Bigbore	21	MFL662B	EVA Infusion Bag 250ml Light Protected Bigbore	21
MF1663	EVA Infusion Bag 500ml with Luerlock	20	MFL663	EVA Infusion Bag 500ml with Luerlock Light Protected	20
MF1663B	EVA Infusion Bag 500ml Bigbore	21	MFL663B	EVA Infusion Bag 500ml Light Protected Bigbore	21
MF1664	EVA Infusion Bag 1000ml with Luerlock	20	MFL664	EVA Infusion Bag 1000ml with Luerlock Light Protected	20
MF1664B	EVA Infusion Bag 1000ml Bigbore	21	MFL664B	EVA Infusion Bag 1000ml Light Protected Bigbore	21
MF1665	EVA Infusion Bag 2000ml with Luerlock	20	MFL665	EVA Infusion Bag 2000ml with Luerlock Light Protected	20
MF1665B	EVA Infusion Bag 2000ml Bigbore	21	MFL665B	EVA Infusion Bag 2000ml Light Protected Bigbore	21
MF1666	EVA Infusion Bag 3000ml with Luerlock	20	MFL666	EVA Infusion Bag 3000ml with Luerlock Light Protected	20
MF1666B	EVA Infusion Bag 3000ml Bigbore	21	MFL666B	EVA Infusion Bag 3000ml Light Protected Bigbore	21
MF1667	EVA Infusion Bag 5000ml with Luerlock	20	MF8881	Infusion Line for Taxol® with 0.2µm Bacteria Filter	24
MF1675	EVA 2-Chamber Bag for TPN (1500 and 500ml)	22			
MF1676	EVA 2-Chamber Bag for TPN (2500ml and 500ml)	22			
MF1821	Bacteria Filter 0.2µm	25			
MF1823	Bacteria Filter 0.2µm, Vented	30			
MF1824	Vented Micro Filter 1.2µm	27			
MF1825	1.2µm Vented Filter	25			
MF1826	Vented Flat Filter 0.2µm (Positive +)	26			
MF1827	Bacteria Filter 0.2µm, Vented Filter (Positive +)	29			
MF1828	Vented Bacteria Filter 0.2µm (+) small	28			
MF1864	EVA Infusion Bag 1000ml with with 3-way Transfer Set	22			
MF1865	EVA Infusion Bag 2000ml with 3-way Transfer Set	22			

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**Fagron Iberica, S.A.U.** · Josep Tapiolas 150
**Fagron SAS** · 19, Bvd Malesherbes · 75008 Paris
**Fagron Nordic A/S** · Kigkurren 8M 2.